

# Installing the Particle Solutions Evaluation License

## Step 1.

Plug the USB stick provided by Brookhaven Instruments into the computer that has Particle Solutions installed.

## Step 2.

Open the file directory of the USB stick. Double click on .txt file shown. This is your evaluation license code.


Name	Date modified	Type	Size
Adobe Reader	11/22/2016 12:32 ...	File folder	
Manuals	9/23/2019 3:28 PM	File folder	
MSDS	12/6/2018 9:03 AM	File folder	
60DayEvaluationLicense.txt	7/16/2019 4:15 PM	Text Document	1 KB
Particle Solutions Software Video .mp4	7/23/2018 3:06 PM	MP4 Video	687,349 KB
PS_3.6.0.7122_Setup.exe	2/3/2020 9:27 AM	Application	40,243 KB

## Step 3.

Select the entire license code string, making sure all characters are selected. You can easily select all of the characters by pressing “Ctrl” and “A” at the same time. Copy the code, either by right-clicking and copying it, or by pressing “Ctrl” and “C” at the same time.

## Step 4.

Run the Particle Solutions Software. Once the software has started, you should see a window similar to the one shown here. Select “Enter License...” which will open a new window.

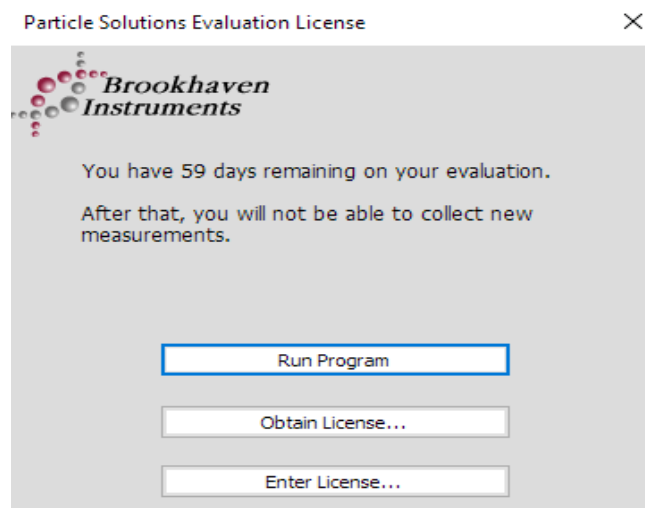

**Step 5.**

Paste your copied license code into the license window, either by right-clicking in the text field and pasting, by using the “Paste From Clipboard” button, or by left-clicking in the text field and pressing “Ctrl” and “V”.

License ✕

**Current License**

Customer Name=Evaluation Customer  
Customer Address=<Evaluation - Unknown>  
Instrument Serial Number = <BIC-Manufacture >  
License Version=3.0


**New License**

lgIAACEORhATPNUBPACZAEN1c3RvbWVyIE5hbWU9RXZhbHV0YXRpb24gQ3VzdG9tZXIjQ3  
VzdG9tZXIqQWRkcmVzcz  
08RXZhbHVhdGlvbiAtIFVua25vd24+I0luc3RydW 1lbnQGU2VyaWFsIE5 1bWJlcj08QkdlLU 1hb  
.....

Customer Name=Evaluation Customer  
Customer Address=<Evaluation - Unknown>  
Instrument Serial Number = <BIC-Manufacture >  
License Version=3.0  
Model=<Unknown>

**Step 6.**

Once you have pasted in your license code string, click on “Apply New License.” You will see a pop-up confirmation window, and the License window will update to say “New license applied.”

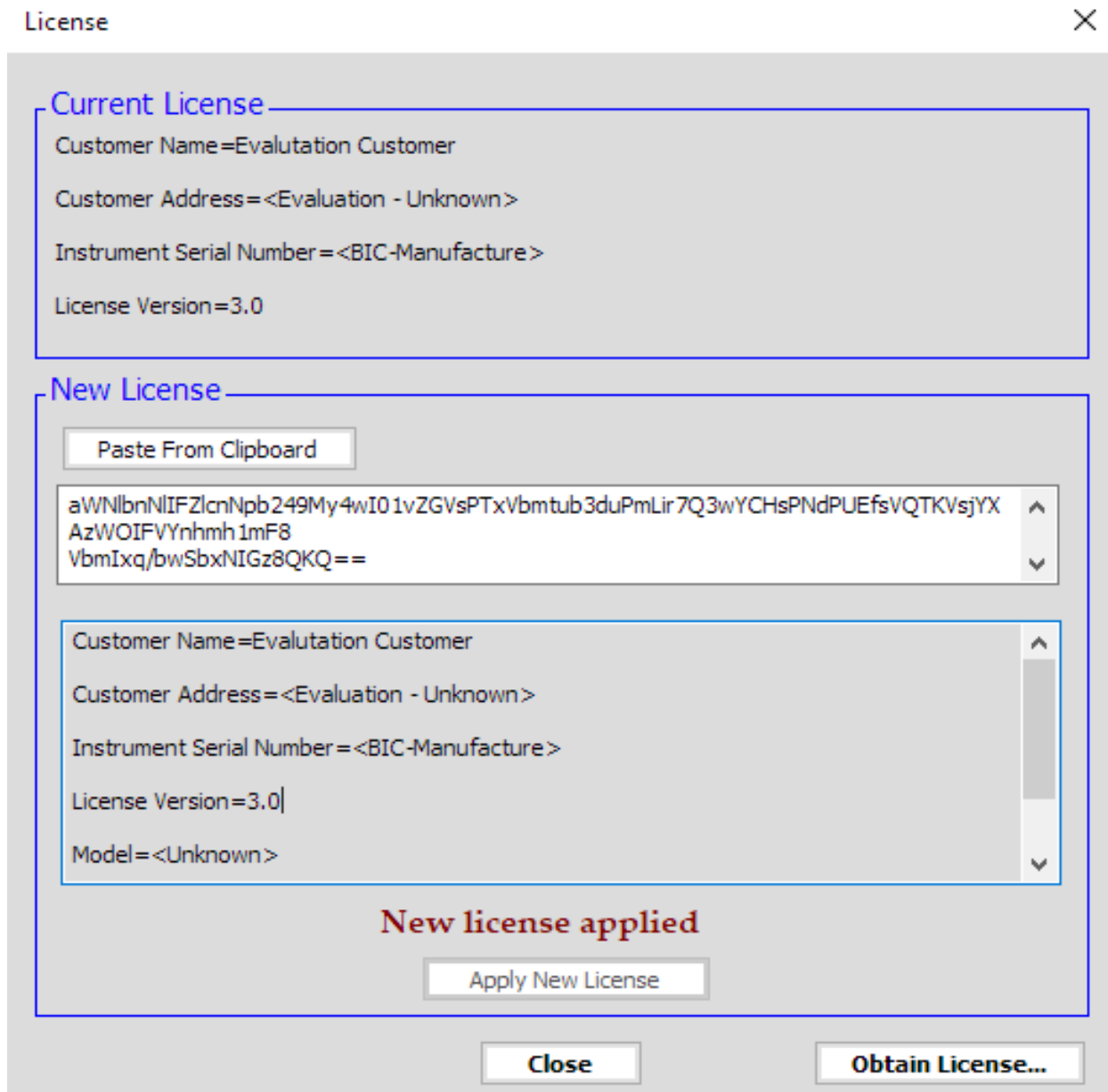

The screenshot shows a window titled "License" with a close button (X) in the top right corner. The window is divided into two main sections: "Current License" and "New License".

**Current License**

- Customer Name=Evaluation Customer
- Customer Address=<Evaluation - Unknown>
- Instrument Serial Number = <BIC-Manufacture >
- License Version=3.0

**New License**

Paste From Clipboard

aWNlbnNlIFZlcnNpb249My4wIO1vZGVsPTxVbmtub3duPmLir7Q3wYCHsPNdPUEfsVQTKVsYX  
AzWOIFVYnhmh1mF8  
VbmIxxq/bwSbxNIGz8QKQ==

Customer Name=Evaluation Customer  
Customer Address=<Evaluation - Unknown>  
Instrument Serial Number = <BIC-Manufacture >  
License Version=3.0|  
Model=<Unknown>

**New license applied**

Apply New License

Close Obtain License...

You may now close the license window and run Particle Solutions.