

Installing the Particle Solutions License

Step 1.


You should receive two attachments with this email. One attachment is this instructional document. The other attachment is a .txt file containing your license code. Open the .txt file to access the license code.

Step 2.

Select the entire license code string, making sure all characters are selected. You can easily select all of the characters by pressing “Ctrl” and “A” at the same time. Copy the code, either by right-clicking and copying it, or by pressing “Ctrl” and “C” at the same time.


Step 3.

Run the Particle Solutions Software. Once the software has started, you should see a window similar to the one shown here. Select “Enter License...” which will open a new window.

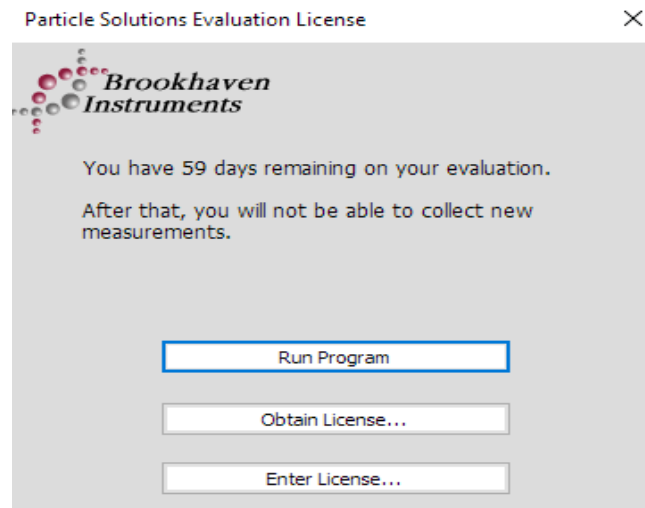

Step 5.

Once you have pasted in your license code string, click on “Apply New License.” You will see a pop-up confirmation window, and the License window will update to say “New license applied.”

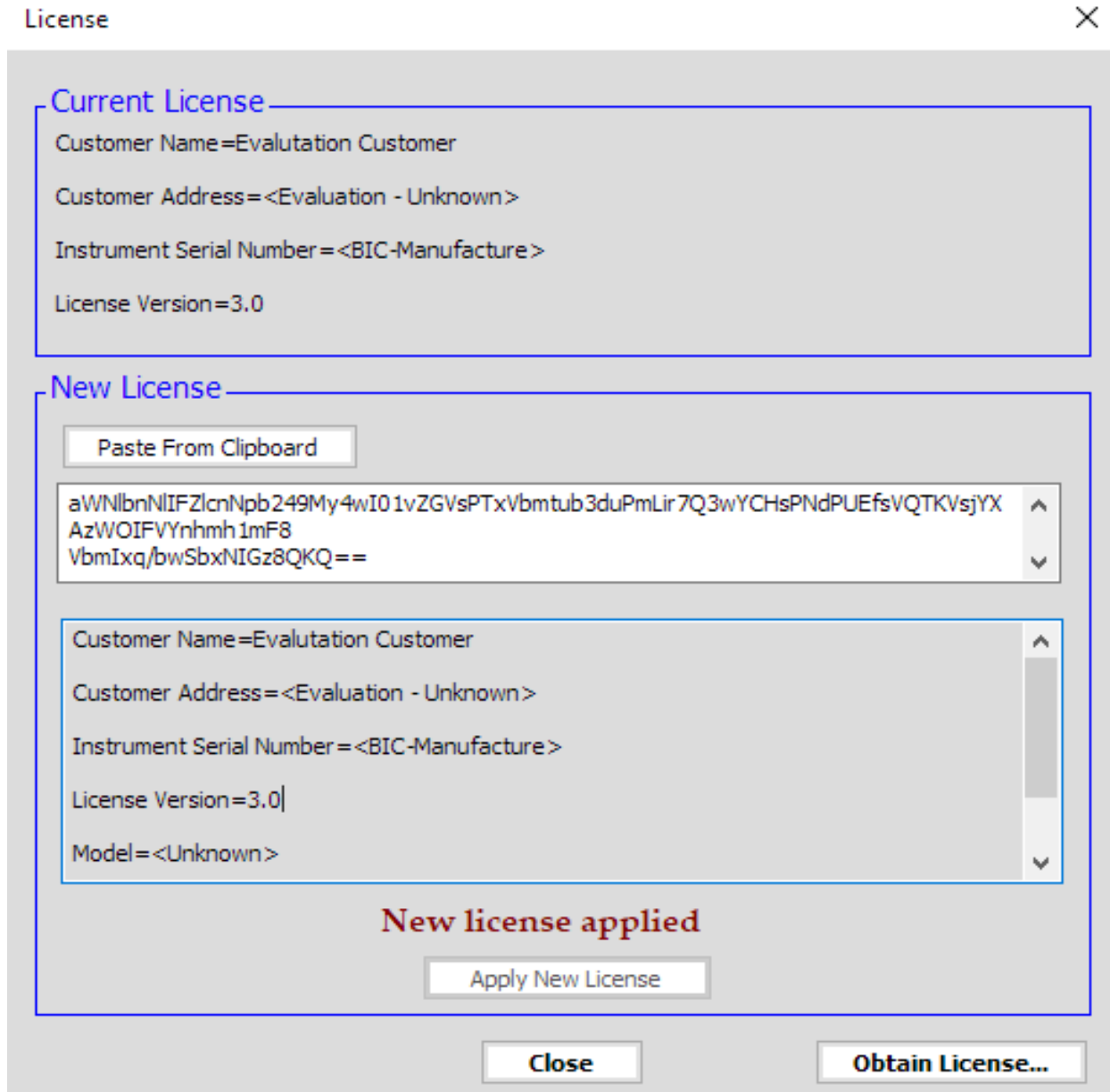

The screenshot shows a window titled "License" with a close button (X) in the top right corner. The window is divided into two main sections: "Current License" and "New License".

Current License:

- Customer Name = Evaluation Customer
- Customer Address = <Evaluation - Unknown >
- Instrument Serial Number = <BIC-Manufacture >
- License Version = 3.0

New License:

- A button labeled "Paste From Clipboard" is located above a text input field.
- The text input field contains the license code: `aWNlbnNlIFZlcnNpb249My4wI01vZGVsPTxVbmtub3duPmLir7Q3wYCHsPNdPUEfsVQTKVsYXAzWOIFVYnhmh1mF8VbmIxcq/bwSbxNIGz8QKQ==`
- Below the input field is a list of license details:
 - Customer Name = Evaluation Customer
 - Customer Address = <Evaluation - Unknown >
 - Instrument Serial Number = <BIC-Manufacture >
 - License Version = 3.0
 - Model = <Unknown >

At the bottom of the "New License" section, the text **New license applied** is displayed in red. Below this text is an "Apply New License" button.

At the bottom of the window, there are two buttons: "Close" and "Obtain License..."

You may now close the license window and run Particle Solutions.